



PFC Sampling Information

ESSENTIALS:

- Wash your hands before opening any of the sampling sets.
- Use the gloves that were provided in the bag.
- Make sure to complete all labeling/paperwork and please be legible.
- Once the appropriate bottles are filled, place them back in the bag that was provided.
- Bottles should be kept on ice until pick up/drop-off.



Drinking Water PFC Set

Method 533

The DW PFC set comes with a total of seven bottles; three 250ml Ammonium Acetate bottles and four unpreserved 250ml Field Blank bottles.

You will fill the three Ammonium Acetate bottles with sampling water.

Then, you will take your unpreserved reagent water and transfer it to the empty Field Blank bottle with the blue label. You will do this again for your second Field Blank. You will then discard the empty bottles from which you poured.

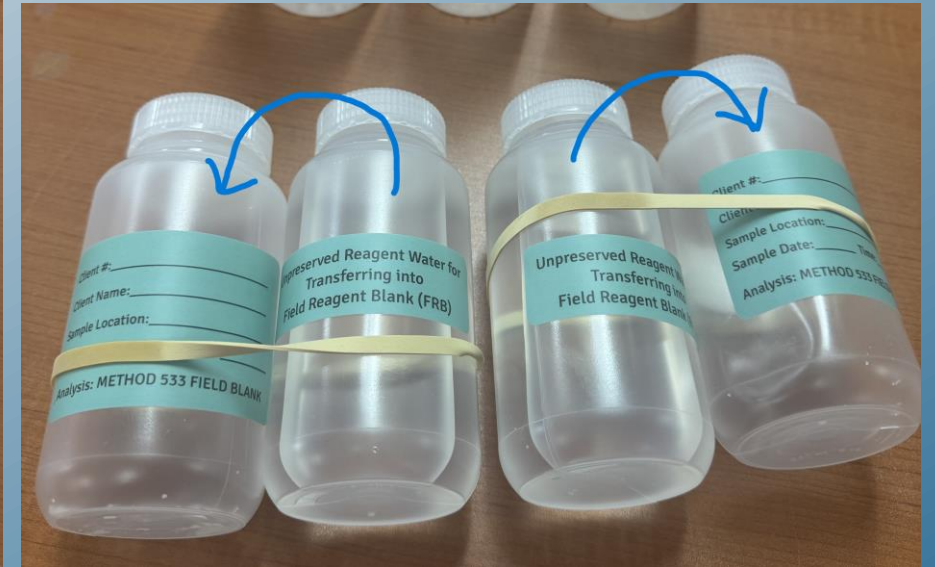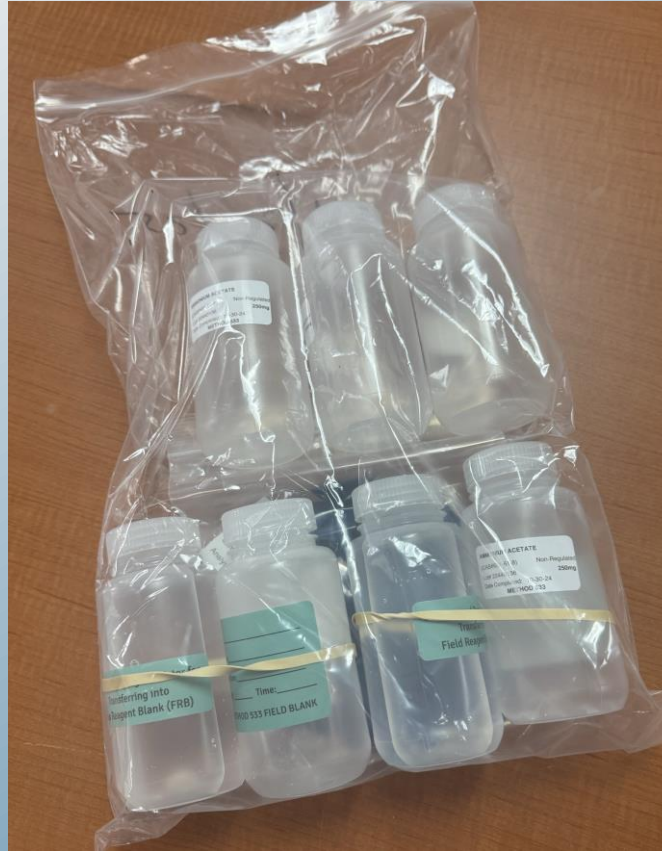

Wastewater PFC Set

Method 1633



The WW PFC set comes with three unpreserved bottles.

All three bottles should be filled to the shoulder for complete sampling.

Sludge PFC Set

Method 1633

The Sludge PFC set comes with two small bottles. Both shall be filled to complete the sampling procedure.

